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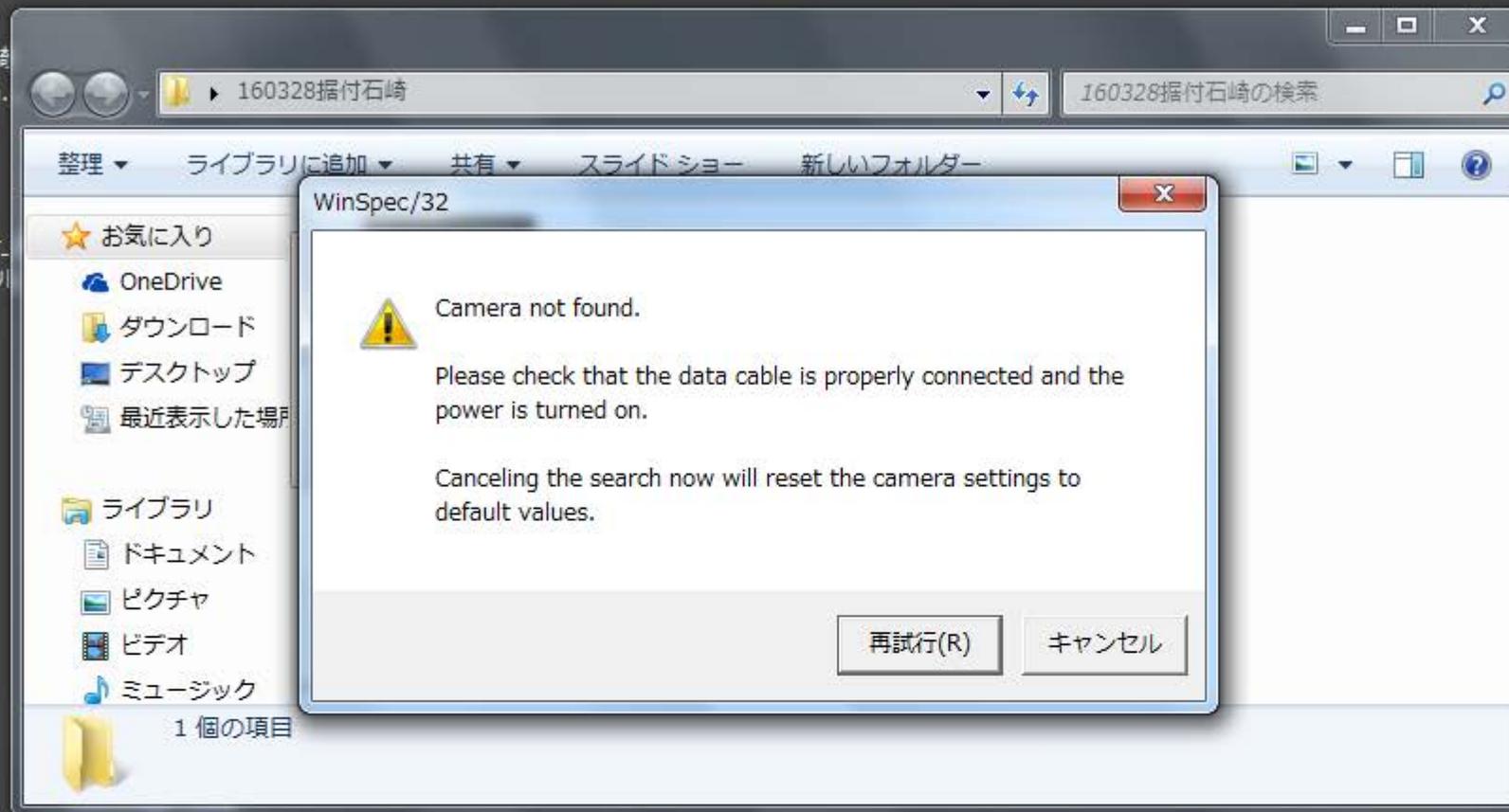
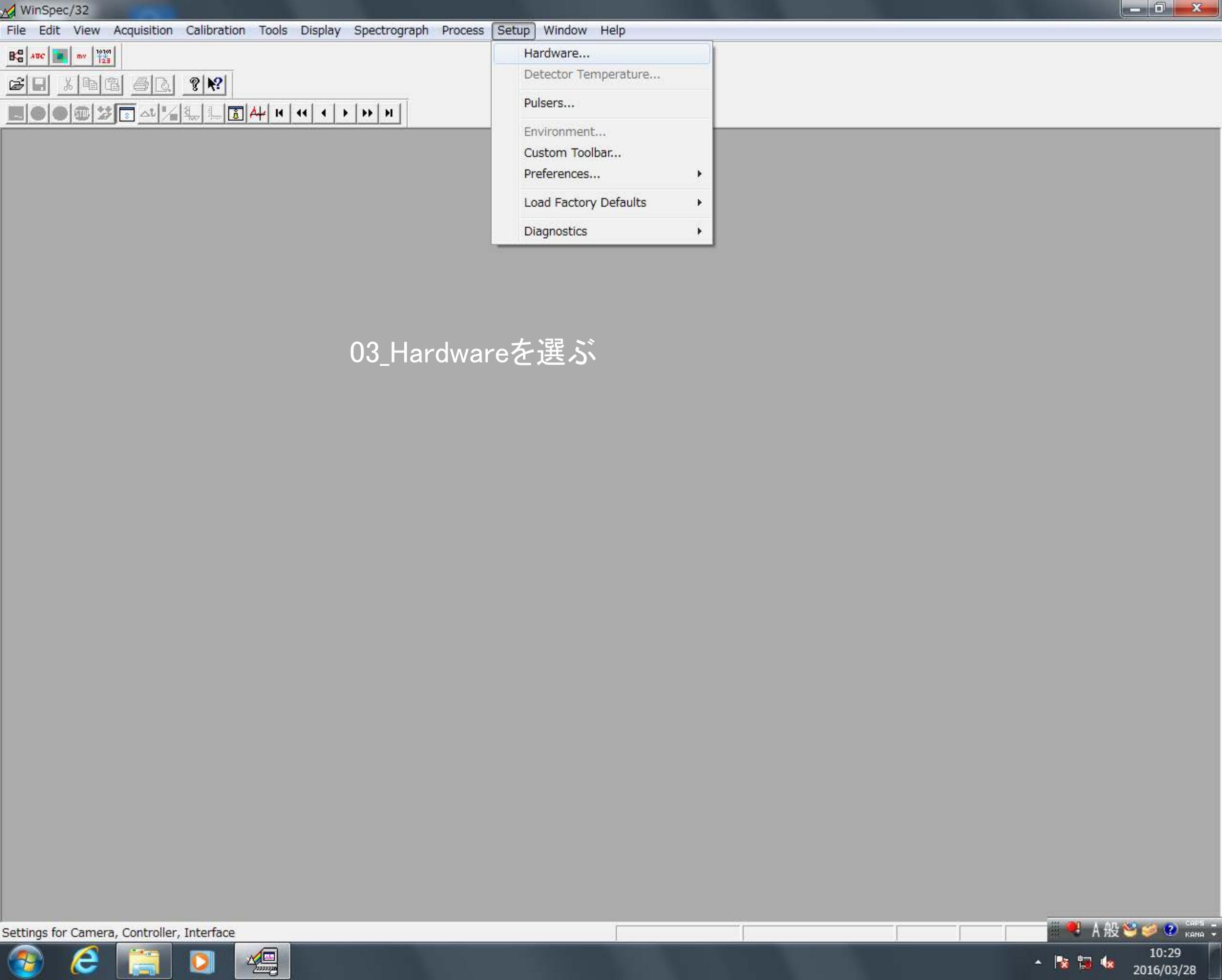
03\_Hardwareを選ぶ

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Camera Detection Wizard - Welcome

This wizard will attempt to detect the camera system you are using. Please make sure your system is connected to the computer and powered on before proceeding.

I will configure the camera system manually

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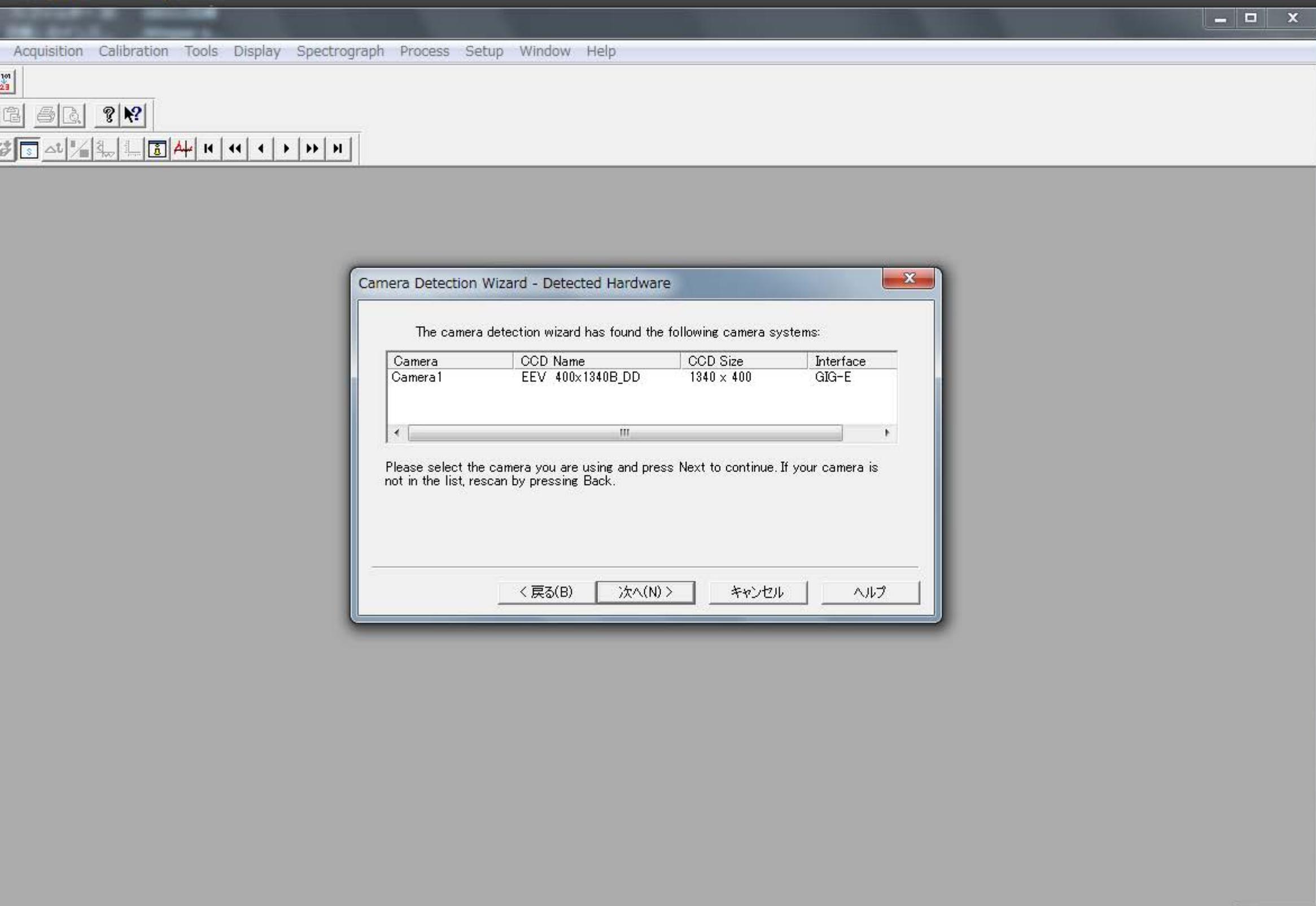
Camera Detection Wizard - Detected Hardware

The camera detection wizard has found the following camera systems:

Camera	CCD Name	CCD Size	Interface
Camera1	EEV 400x1340B_DD	1340 x 400	GIG-E

Please select the camera you are using and press Next to continue. If your camera is not in the list, rescan by pressing Back.

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System tray area showing system icons (network, volume, power), user name (A 般), and date/time (10:36, 2016/03/28).

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Acquisition Calibration Tools Display Spectrograph Process Setup Window Help

101  
23

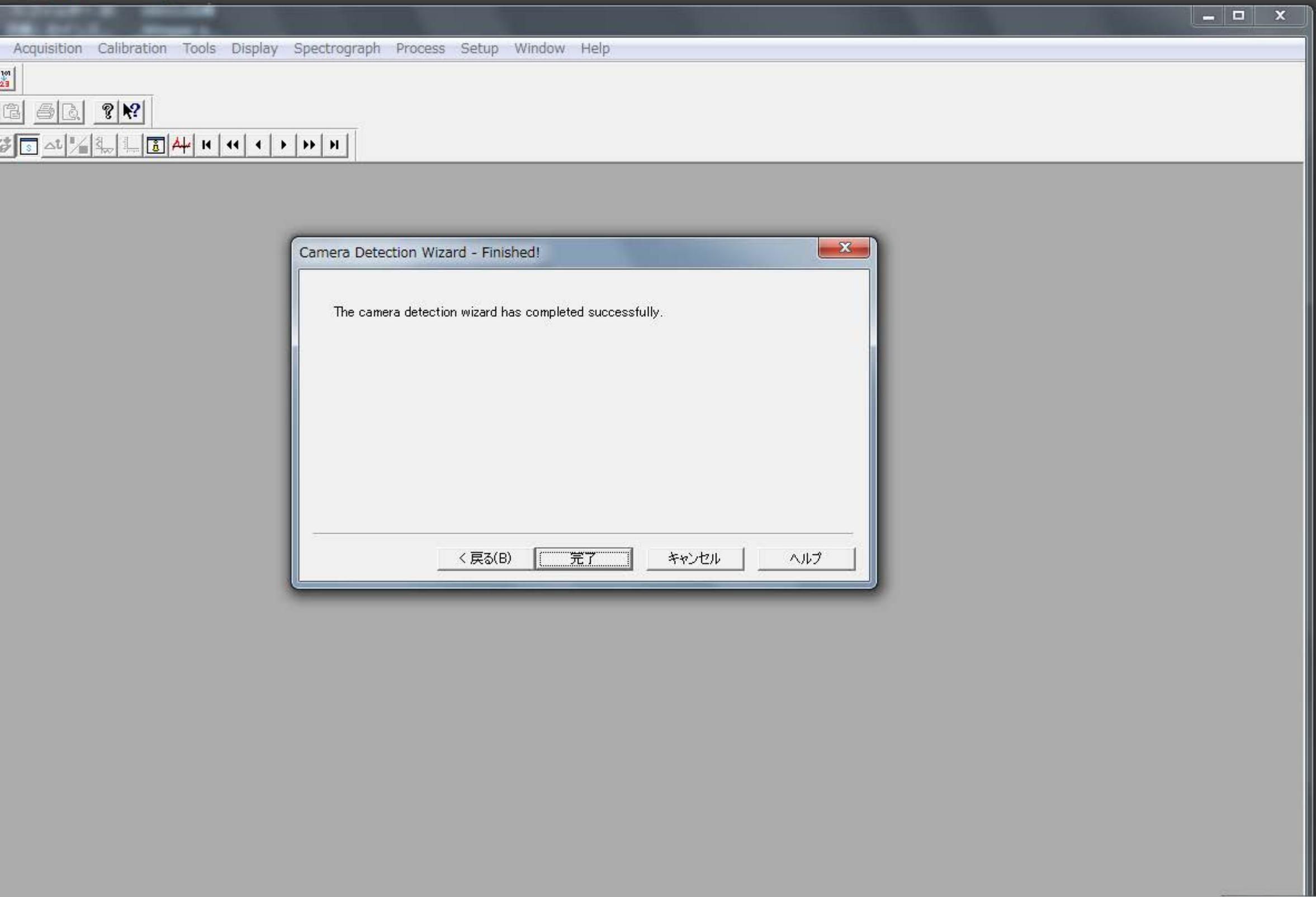
Camera Detection Wizard - Test Image

Would you like to aquire a test image?

Yes  No

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WinSpec/32

File Edit View Acquisition Calibration Tools Display Spectrograph Process Setup Window Help

ABC mv 1010 123

File Edit View Acquisition Calibration Tools Display Spectrograph Process Setup Window Help

File Edit View Acquisition Calibration Tools Display Spectrograph Process Setup Window Help

Move ...

Define ...

Calibrate...

A 般 CAPS KANA

10:40  
2016/03/28

WinSpec/32

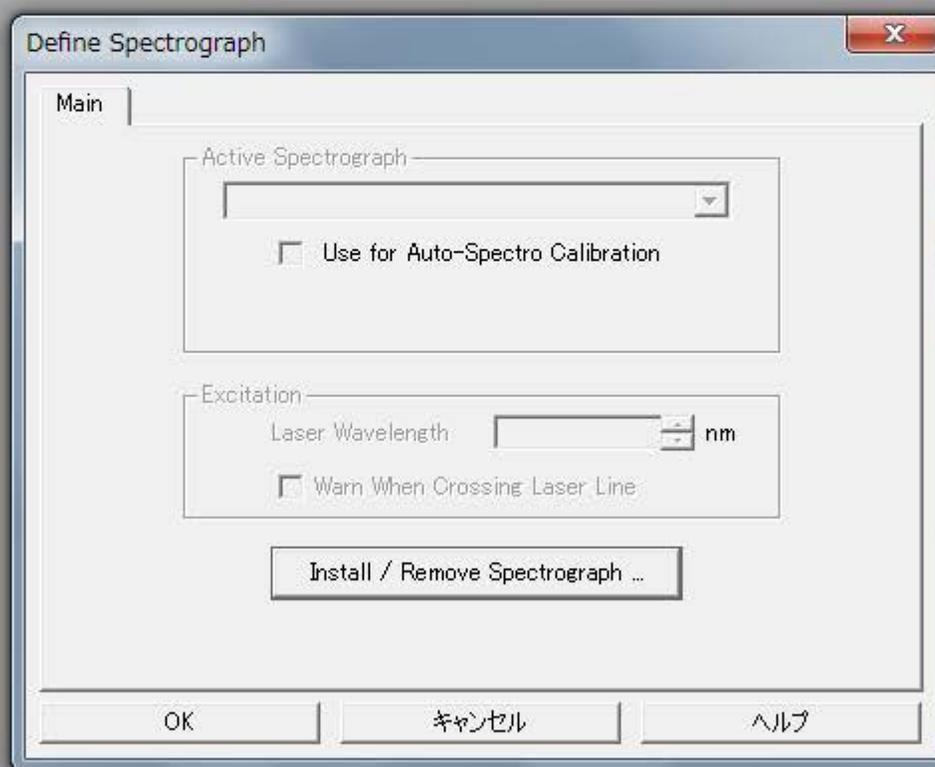
File Edit View Acquisition Calibration Tools Display Spectrograph Process Setup Window Help

ABC mv 1010 123

File Edit View Acquisition Calibration Tools Display Spectrograph Process Setup Window Help


File Edit View Acquisition Calibration Tools Display Spectrograph Process Setup Window Help

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
CAPS KANA

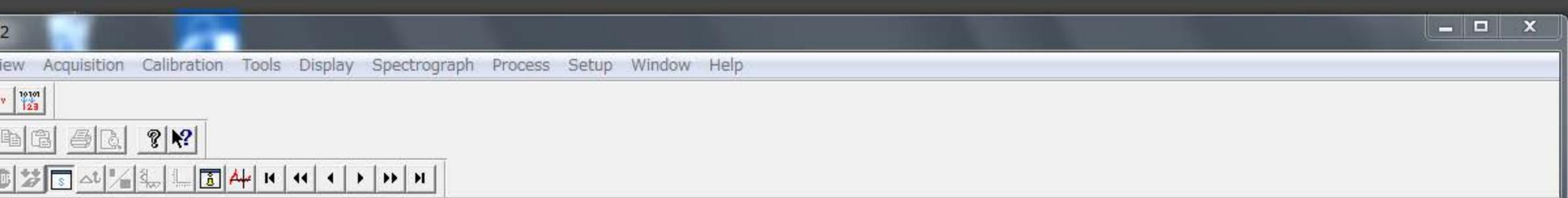10:41  
2016/03/28

## 10\_Acton SP500iを選ぶ

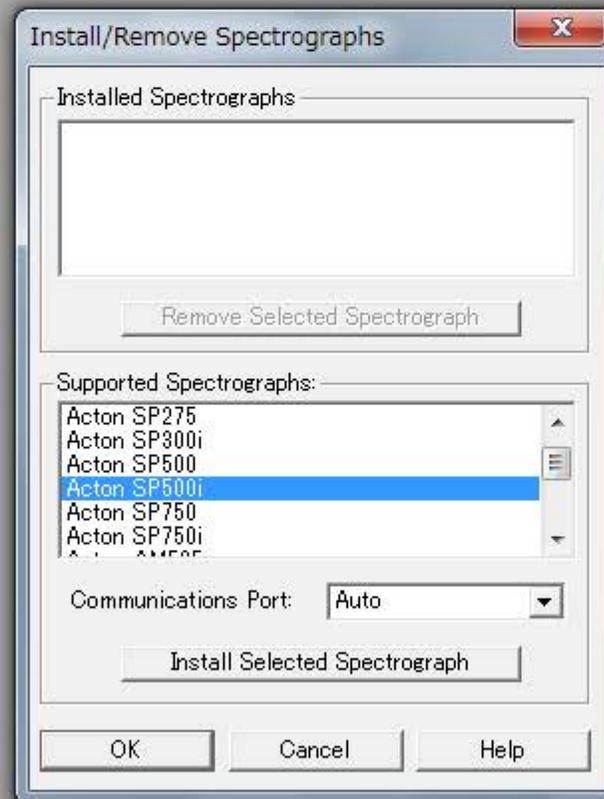WinSpec/32

File Edit View Acquisition Calibration Tools Display Spectrograph Process Setup Window Help

ATC mv 1010 123

File Edit View Acquisition Calibration Tools Display Spectrograph Process Setup Window Help

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